


XRF (Malvern Panalytical, Epsilon 3)

Cautions on the usage 利用上の注意点

During measurement, your sample cup is set on the measurement chamber as shown in the below image.

測定中は、下図のように試料カップが測定部にセットされます。

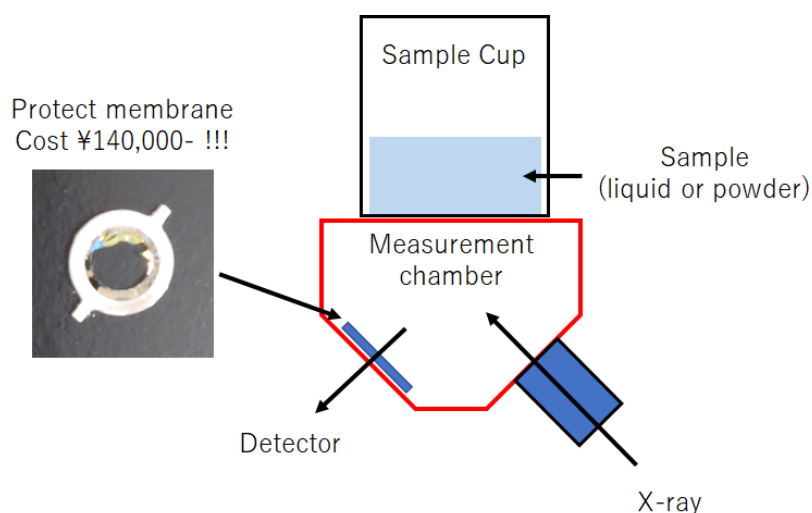

If the sample falls into the chamber, the machine got error and the protect membrane may be damaged. So, please endure the following.

1. Before measurement, please carefully check that the sample cup has no leakage, and the bottom part is clean.
2. When you find the cup has been damaged after measurement, please immediately let me know (ex. 4636, toruwada@jaist.ac.jp).

試料がチャンバー内に落下すると、機械のエラーや保護膜の破損の原因となります。そのため、以下の点にご注意ください。

1. 測定前に、試料カップに液漏れがないこと、下面が汚れていないことをよく確認してください。
2. 測定後、カップが破損しているのを発見した場合、直ちにご連絡ください（内線：4636、toruwada@jaist.ac.jp）。